



Laser Eyewear Cleaning Procedure

Laser eyewear should be cleaned with a mild detergent soap and plenty of lukewarm water. Attention should be paid to the parts of the eyewear that directly touch the skin (e.g., bridge of nose, ears) and care should be taken when cleaning and drying the lens to avoid scratches, especially with glass eyewear with dielectric coatings.

AVOID harsh chemicals and other cleaning products as they may damage the lens. If your lab does not have an internal sink, please clean the eyewear in a restroom.

Email radiation_safety@harvard.edu to send comments and suggestions to the EH&S Radiation Safety Office.